
INSTRUCTIONS FOR ASSEMBLY

MODEL C-10 SERIES CHROMATO-VUE® CABINET

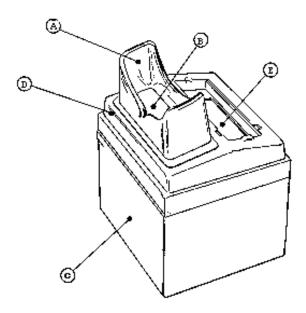


- 1. Lift flexible eyepiece (A) and cover assembly (D) out through top of cabinet.
- Slip lamp well (E) portion of cover assembly over rim at back of cabinet until flush with top. Apply enough pressure toward front portion to snap over rim. Ensure that cover assembly is flush with top on all sides.
- 3. Slip flexible eyepiece over rim above contrast control filter (B) with notched portion facing front of cabinet.
- 4. Place the Mineralight® lamp into the molded lamp well with power cord placed to the left. The Chromato-Vue cabinet is now ready to use.
- 5. Lift light curtain (C) when placing or positioning materials; leave in closed position when in use to block external light from entering the cabinet.



INSTRUCTIONS FOR ASSEMBLY

MODEL C-10 SERIES CHROMATO-VUE® CABINET

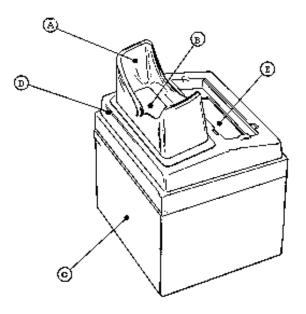

- 1. Lift flexible eyepiece (A) and cover assembly (D) out through top of cabinet.
- 2. Slip lamp well (E) portion of cover assembly over rim at back of cabinet until flush with top. Apply enough pressure toward front portion to snap over rim. Ensure that cover assembly is flush with top on all sides.
- 3. Slip flexible eyepiece over rim above contrast control filter (B) with notched portion facing front of cabinet.
- 4. Place the Mineralight® lamp into the molded lamp well with power cord placed to the left. The Chromato-Vue cabinet is now ready to use.
- 5. Lift light curtain (C) when placing or positioning materials; leave in closed position when in use to block external light from entering the cabinet.



CHROMATO-VUE CABINET C-10P

bracket is to hold the portable lamp securely in place. addition of a lamp support bracket on the C-10P. The purpose of the The difference between the C-10, C-10E and C-10P Cabinets is the

C-10P is as follows: numbers 1 through 3 on the previous page. Instruction number 4 for the To assemble the C-10P, follow the C-10 Cabinet Series instructions

block out any outside light. materials; leave the curtain in the closed position when in use to ready to use. Lift light curtain (C) when placing or positioning firmly holds the lamp in place. The Chromato-Vue cabinet is now 4. Place the portable lamp in the lamp support bracket. This bracket

Replacement Parts:

10-6100-86 Contrast Control Filter: 23-0036-01

Eyepiece:

(606) 646-3197 / Fax: (909) 946-3597 2066 W. 11th Street Upland, CA 91786 / Tel: (800) 452-6788 Corporate Headquarters: UVP, LLC

Eax: +44(0)1223 420156 Cambridge CB4 1TG UK / Tel: +44(0)1223 420022 Unit 1, Trinity Hall Farm Estate, Nuffield Road, European Operations: Ultra-Violet Products Ltd

For additional information on UVP's products, go to www.UVP.com

1-0008-01 Rev J

Mineralight and Chromato-Vue are registered trademarks of UVP, LLC

CHROMATO-VUE CABINET C-10P

bracket is to hold the portable lamp securely in place. addition of a lamp support bracket on the C-10P. The purpose of the The difference between the C-10, C-10E and C-10P Cabinets is the

C-10P is as follows: numbers 1 through 3 on the previous page. Instruction number 4 for the To assemble the C-10P, follow the C-10 Cabinet Series instructions

block out any outside light. materials; leave the curtain in the closed position when in use to ready to use. Lift light curtain (C) when placing or positioning firmly holds the lamp in place. The Chromato-Vue cabinet is now Place the portable lamp in the lamp support bracket. This bracket

Eyepiece: 23-0036-01 Replacement Parts:

38-0013-01 Contrast Control Filter:



7638-349 (609) 346-349 (609) 2066 W. 11th Street Upland, CA 91786 / Tel: (800) 452-6788 Corporate Headquarters: UVP, LLC

Fax: +44(0)1223 420156 Cambridge CB4 1TG UK / Tel: +44(0)1223 420022 Unit 1, Trinity Hall Farm Estate, Nuffield Road, European Operations: Ultra-Violet Products Ltd

For additional information on UVP's products, go to www.UVP.com

しv9月 f0-8000-f8

Mineralight and Chromato-Vue are registered trademarks of UVP, LLC