


Ultraviolet Blocking Eyewear

UVC-503 Goggles, UVC-303 Spectacles and UVC-803 Faceshield

Eyewear block ultraviolet radiation while providing optimum contrast in viewing and lessened eye fatigue. The UVC-503 Goggles have air vents that are located on each side of the goggles to allow air flow while blocking the ultraviolet radiation. The UVC-303 Spectacles are lightweight and fit comfortably on the face, even over prescription glasses. Side arm panels give additional eye protection. The UVC-803 Faceshield covers the neck and face area.

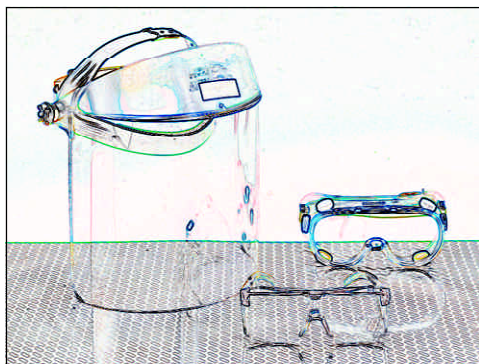


Important: The eyewear are intended to protect the eyes from harmful ultraviolet radiation emitted from UV lamps. The eyewear is designed for eye protection from lamps emitting wavelengths of 254nm to 365nm (nanometers). These Ultraviolet Blocking Eyewear also reduce longwave ultraviolet “blue haze” for optimum viewing.

For exposure periods other than eight hours or for ultraviolet wavelength emissions other than 254nm to 365nm, consult Document HSM 73-11009 “Criteria for a Recommended Standard ... Occupational Exposure to Ultraviolet Radiation published by the US Department of Health, Education and Welfare.”

Caution: Although the Goggles and Spectacles are intended to protect the eyes and the Faceshield to protect the face and neck, care must be taken to protect other parts of the body from severe skin burn associated with short-wave UV wavelengths. Avoid shining any ultraviolet lamp on reflective surfaces. Wear protective clothing and skin cream.

Eyewear meet performance standards of ANSI Z87.1-1989. Spectacles and Goggles lenses are impact resistant, but not unbreakable. Clean and inspect frequently. Pitted or scratched eyewear reduce protection — replace immediately. Clean all eyewear surfaces with a mild soap solution. Do not use solvent on lenses. Carefully rinse off all soap. Air dry or pat dry with a clean, soft cloth or tissue.

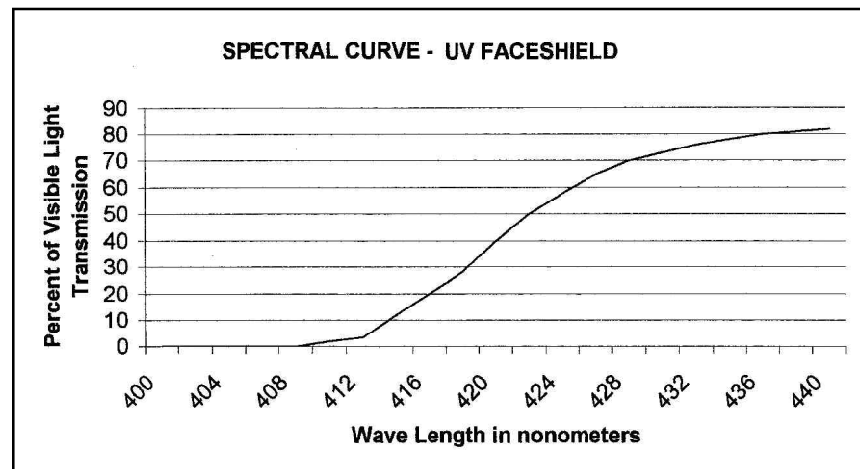

Ultraviolet Blocking Eyewear

Model	Description	Parts Numbers
UVC-303	Spectacles	98-0002-01
UVC-503	Goggles	98-0002-02
UVC-803	Faceshield	98-0002-04
	Replacement lens for Goggles	98-0002-03
	Shield Replacement	38-0151-01

Product Warranty. UVP’s products are guaranteed to be free of defects in materials, workmanship and manufacture for one (1) years from the date of purchase. If equipment failure or malfunction occurs during the warranty period, UVP shall examine the inoperative equipment and have the option of repairing or replacing any part(s) which, in the judgment of UVP, were originally defective or became so under conditions of normal usage and service.



Web Site: www.uvp.com

Corporate Headquarters: UVP, LLC

2066 W. 11th Street, Upland, CA 91786 | Tel: (800)452-6788
or (909)946-3197 | Fax: (909)946-3597

European Operations: Ultra-Violet Products Ltd.

Unit 1, Trinity Hall Farm Estate, Nuffield Road, Cambridge CB4 1TG UK
Tel: +44(0)1223-420022 | Fax: +44(0)1223-420561